



Designed to help you choose
the right membrane filtration unit.

Product Guide

AN	DS	TB	PG	SG
----	----	----	----	----

Drug Compatibility Guide for Disposable Syringe Filters

MILLIPORE



ADVANCING LIFE SCIENCE TOGETHER™
Research. Development. Production.

Millipore and Millex are registered trademarks of Millipore Corporation.
ISO is a registered trademark of the International Standards Organization.
Life Sciences and Advancing Life Science Together are trademarks of Millipore Corporation.
LIFE No. P55024EN00 12/07 Printed in U.S.A. 07-396 © 2007 Millipore Corporation,
Billerica, Massachusetts 01821 U.S.A. All Rights Reserved.

www.millipore.com/offices



Choosing the right membrane filtration unit.

Millipore's products have become the standard in the medical industry due to their consistent high quality and reliable performance. Many hospitals, pharmacies, clinical laboratories, and home health care facilities around the world have relied on Millipore for over 50 years.

The Millex® Filter Unit, designed for small-volume sterilization and clarification of pharmacy solutions, is our most popular filtration product. Like all Millipore filters, Millex units are built with quality in mind. Raw materials are carefully inspected before being used and we exercise stringent control over our manufacturing process. Our production facility complies with FDA Good Manufacturing Practices (cGMP) and has earned ISO® 13485 certification specific for medical devices. Numerous quality checks are done to make sure that our products will work as specified time after time. Our attention to detail throughout the whole production process is what separates Millipore filters from the rest.



Backed by Knowledgeable Technical Support

Not only do we supply high quality products, like Millex filter units, but we also have a staff of technical representatives available to answer any questions you may have about your particular filtration application. Just pick up the phone and call **1-800-MILLIPORE** (1-800-645-5476). Our representatives can help you select the product you need or assist you in resolving a problem. For information outside of North America, please visit www.millipore.com/offices.

Millipore is dedicated in developing products and providing services that improve productivity in biopharmaceutical manufacturing and in clinical, analytical and research laboratories. Visit www.millipore.com to see all of our comprehensive products and services.

Filter Compatibility Chart

Drug Name	Millex GV	Millex GS	Millex OR	Millex FG	Millex HA	Millex HV	Millex SV	Reference
5-Fluorouracil			○	○				30, 31
[18F] Fluoromisonidazole, Misiomidazole	○							5
Acebutolol		○						12
Acetylsalicylic acid		○						12
Alpha1-Proteinase Inhibitor (Human)							○	28
Alprenolol		○						12
Amiloride		○						12
Amphotericin B for Injection USP							○	14, 21, 24, 32, 37
Atenolol		○						12
Azathioprine					○	○		35
Azodicarbonamide		○						6
Bleomycin Sulfate			○	○				30, 31
Caffeine		○						12
Centirizine					○	○		34
Chlorothiazide		○						12
Chloramphenicol		○						12
Cimetidine		○						12
Ciprofloxacin		○						29
Cisplatin, Cisplatin Injection			○	○				30, 31
Cyclosporine A	○							10
Cytarabine			○	○				30
Daunorubicin			○	○				2, 30, 31
DE-310		○						27
Diclofenac						○		1
Enalaprol		○						12
Ethionamide				○				17
Factor IX Complex Heat-Treated							○	28
Gatifloxacin					○	○		22
Hydrochlorothiazide		○						12
Ibuprofen					○	○		1
Isoniazide				○				17
isonicotinic acid				○				17
Ketamine		○						26
Las 35917						○		15
Levofloxacin					○	○		22
Lomefloxacin					○	○		22
Methyl Gag; NSC-32946			○					30, 31
Metoprolol		○						12
Mitomycin			○	○				30, 31
Morphazinamide				○				17
Nadolol		○						12
Nicotinic acid				○				17
Paclitaxel	○							7
p-Aminobenzoic acid (PABA)						○		1
p-aminosalicylic acid				○				17
Pefloxacin					○	○		22
Pentoxifylline (PTX)	○							23
Phenytoin							○	33
Pyrazinamide				○				17
Pyrimethamine					○	○		20
Ranitidine		○						12
Rifampicin					○	○		16
Sabeluzole						○		11
Streptokinase							○	24
Sulfadozine						○		20
Sulphasalazine		○						12
Sulpride		○						12
Terbutaline		○						12
Thiotepa Parenteral Sterile			○	○				30, 31
Timolol		○						12
Tobramycin Vincristine Sulfate			○	○				4, 8, 9, 13, 18, 19, 30, 31, 36
Tranexamic acid		○						12
Triamcinolone Acetonide		○						25
Triazine; NSC-139105			○	○				30
Tropicamide					○			3
Vinblastine Sulfate				○				30, 31

References

- Santini, A. O., et al. (2004). p-Aminobenzoate ion determination in pharmaceutical formulations by using a potentiometric sensor immobilized in a graphite matrix. *Talanta* **63**: 833-838.
- Bosanquet, A.G. (1985). Stability of solutions of antineoplastic agents during preparation and storage for *in vitro* assays. General considerations, the nitrosoureas and alkylating agents. *Cancer Chemother. Pharmacol.* **14**: 83-95.
- Cappello, B., et al. (2001). Solubilization of tropicamide by hydroxypropyl- β -cyclodextrin and water-soluble polymers: *in vitro/in vivo* studies. *Int. J. Pharm.* **213**: 75-81.
- Butler, L.D., et al. (1980). Effect of inline filtration on the potency of low-dose drugs. *Am. J. Hosp. Pharm.* **37**: 935-941.
- Chang, C.W., et al. (2007). A robotic synthesis of [18F]fluoromisonidazole ([18F]FMISO), *Applied Radiation and Isotopes*, In Press, Accepted Manuscript, Available online 6 February.
- Fagny, C., et al. (2002). Ribonucleotide reductase and thymidine phosphorylation: two potential targets of azodicarbonamide. *Biochem. Pharmacol.* **64**: 451-456.
- Kiesewetter, et al. (2003). Fluoro-, bromo-, and iodopaclitaxel derivatives: synthesis and biological evaluation. *Nuclear Medicine and Biology* **30**: 11-24.
- Elenbaas, J.K., et al. (1985). Effect of inline filtration on tobramycin delivery. *Drug Intell. Clin. Pharm.* **19**: 122-125.
- Ennis, C.E., et al. (1983). *In vitro* study of inline filtration of medications commonly administered to pediatric cancer patients. *J. Parenter. Enter. Nutr.* **7**: 156-158.
- Lallemand, et al. (2005) A novel water-soluble cyclosporine A prodrug: Ocular tolerance and *in vivo* kinetics. *Int. J. Pharm.* **295**: 7-14.
- Verreck, G., et al. (2005). Preparation and physicochemical characterization of biodegradable nerve guides containing the nerve growth agent sabeluzole. *Biomaterials* **26**: 1307-1315.
- Flaten, G.E., et al. (2006). Drug permeability across a phospholipid vesicle based barrier: A novel approach for studying passive diffusion. *European Journal of Pharmaceutical Sciences* **27**: 80-90.
- Grant, A. M. (1987). Personal communication.
- Huber, R.C. and Riffkin, C. (1975). Inline final filters for removing particles from amphotericin B infusions. *Am. J. Hosp. Pharm.* **32**: 173-176.
- Toro, I. et al. (2004). Development and validation of a capillary electrophoresis method with ultraviolet detection for the determination of the related substances in a pharmaceutical compound. *J. Chromatogr. A* **1043**: 303-315.
- Calleja, I., et al. (2004). High-performance liquid-chromatographic determination of rifampicin in plasma and tissues. *J. Chromatogr.* **1031**: 289-294
- Gennaro, M.C., et al. (2001). Ion interaction reagent reversed-phase high-performance liquid chromatography determination of anti-tuberculosis drugs and metabolites in biological fluids. *J. Chromatogr. B* **754**: 477-486.
- Kanke, M., et al. (1983). Binding of selected drugs to a "treated" inline filter. *Am. J. Hosp. Pharm.* **40**: 1323-1328.
- Khue, N.V., and Jung, L. (1985). Study of the retention of child-dose drugs on cellulose ester membranes during inline intravenous filtration. *S-T-P-Pharma* **1**: 201-207.
- Boca, M.B., et al. (2005). A validated HPLC method for determining residues of a dual active ingredient anti-malarial drug on manufacturing equipment surfaces. *J. Pharm. Biomed. Anal.* **37**: 461-468.
- Maddus, M.S., and Barriere, S.L. (1980). A review of complications of amphotericin B therapy: recommendations for prevention and management. *Drug Intell. Clin. Pharm.* **14**: 177-181.
- Ines, M., et al. (2006). Quantitative determination of gatifloxacin, levofloxacin, lomefloxacin and pefloxacin fluoroquinolonic antibiotics in pharmaceutical preparations by high-performance liquid chromatography. *J. Pharm. Biomed. Anal.* **40**: 179-184.
- Fantin, M., et al. (2006). Pentoxifylline and its major oxidative metabolites exhibit different pharmacological properties. *European Journal of Pharmacology* **535**: 301-309.
- McEvoy, G.K. (Ed.). (1991). American hospital formulary service drug information 91. *American Society of Hospital Pharmacists*.
- Ober, M.D. et al. (2006). Measurement of the Actual Dose of Triamcinolone Acetonide Delivered by Common Techniques of Intravitreal Injection. *American Journal of Ophthalmology* **142**: 597-600.e1.
- Lofwall, M.R., et al. (2006). Cognitive and Subjective Acute Dose Effects of Intramuscular Ketamine in Healthy Adults. *Experimental and Clinical Psychopharmacology* **14**: 439-449.
- Kato, M., et al. (2005). Examination of meningocele induced by the antitumor agent DE-310 in rat fetuses. *Reproductive Toxicology* **20**: 495-502. 28. Miles Pharm.
- Langer, O., et al. (2003). Synthesis of fluorine-18-labeled ciprofloxacin for PET studies in humans. *Nuclear Medicine and Biology* **30**: 285-291.
- Pavlik, E.J., et al. (198e). Properties of anticancer agents relevant to *in vitro* determinations of human tumor cell sensitivity. *Cancer Chemother. Pharmacol.* **11**: 8-15.
- Pavlik, E.J., et al. (1982). Sensitivity to anticancer agents *in vitro*: standardizing the cytotoxic response and characterizing the sensitivities of a reference cell line. *Gynecol. Oncol.* **14**: 243-261.
- Oradell, N.J. (1991). *Physicians' Desk Reference*. 45th edition.
- Rusmin, S., et al. (1977). Effect of inline filtration on the potency of drugs administered intravenously. *Am. J. Hosp. Pharm.* **34**: 1071-1074.
- Rudaz, S., et al. (2003). Veuthey, Development and validation of a heart-cutting liquid chromatography-mass spectrometry method for the determination of process-related substances in cetirizine tablets. *Analytica Chimica Acta* **492**: 271-282.
- Fazio, T.T., et al. (2007). Quantitative determination and sampling of azathioprine residues for cleaning validation in production area. *J. Pharm. Biomed. Anal.* **43**: 1495-1498.
- Thompson, D.F., et al. (1984). Effect of inline filtration on pediatric doses of gentamicin and tobramycin. *Infusion* **8**: 31-32.
- Tipple, M., et al. (1977). Availability of active amphotericin B after filtration through membrane filters. *Am. Rev. Resp. Dis* **115**: 879-881.

